

LEXT OLS4100
Offline software

Installation Manual

Installation of OLYMPUS LEXT OLS4100 (Offline Connection)

1.1. Notes

- Offline Connection must be installed according to steps of [2.2] Installation of OLYMPUS LEXT OLS4100.
- During installation, if the following Windows Update screen appears, please select [Postpone].

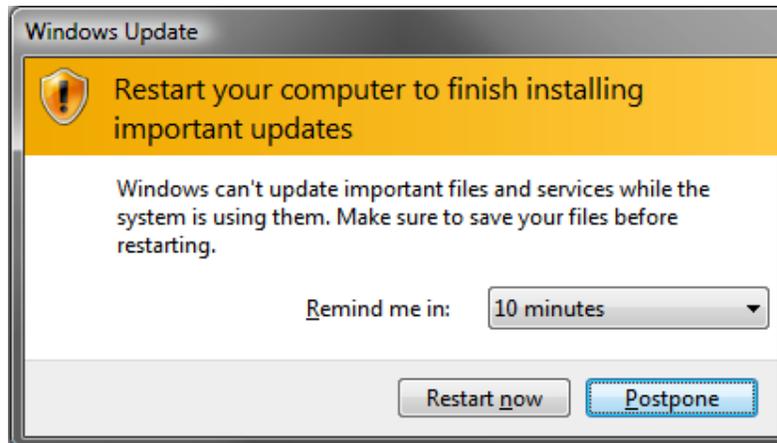


Fig.1 Windows Update

(If this screen appears, select [Postpone].)

- During installation, if Program Compatibility Assistant screen appears shown below, please select [This program was installed correctly].
(Even if the following screen is displayed, the installation is correctly done.)

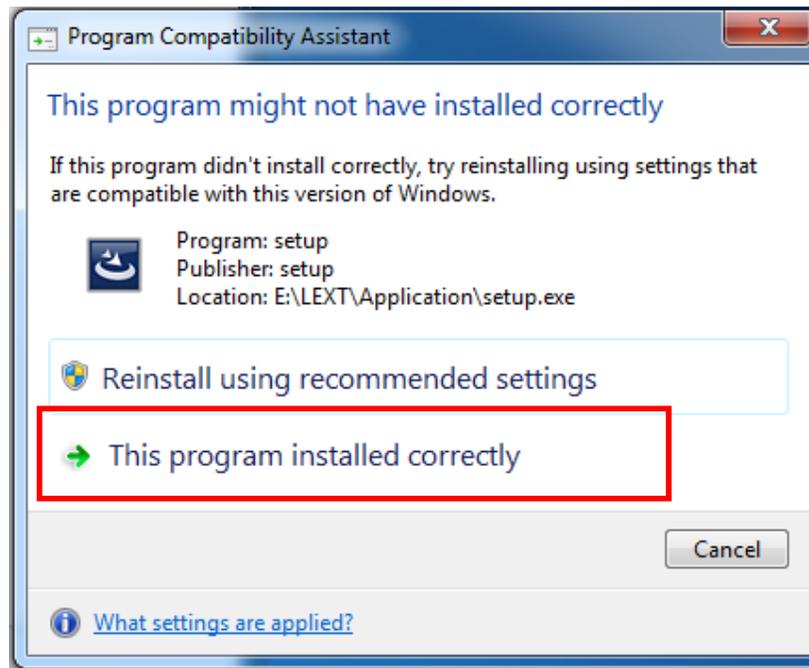


Fig.2 Program Compatibility Assistant screen

(If this screen appears, select **[This program was installed correctly.]**)

- Please remove the exclusive USB dongle key prior to installation.

1.2. Installation of OLYMPUS LEXT OLS4100

*** Please check that the exclusive USB dongle key is not connected.**

If the exclusive USB dongle key is connected, remove it before starting installation.

Perform steps described below.

- ① Insert the installation medium into Drive.
- ② When PC is 32 bit, right-click setup.exe (or setup) in the folder of "LEXT" in a "32bit" folder, and left-click [Run as Administrator (A)] in the menu displayed.
When PC is 64 bit, right-click setup.exe (or setup) in the folder of "LEXT" in a "64bit" folder, and left-click [Run as Administrator (A)] in the menu displayed.
- ③ Operate according to the figures shown below.

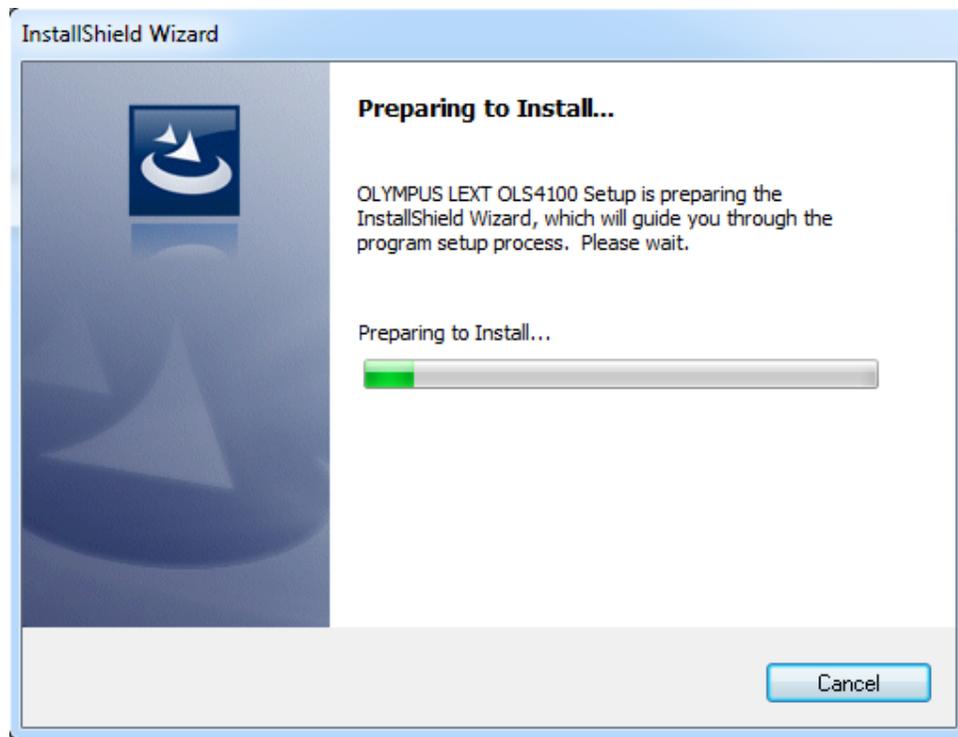


Fig.3 Preparation to install OLYMPUS LEXT OLS4100



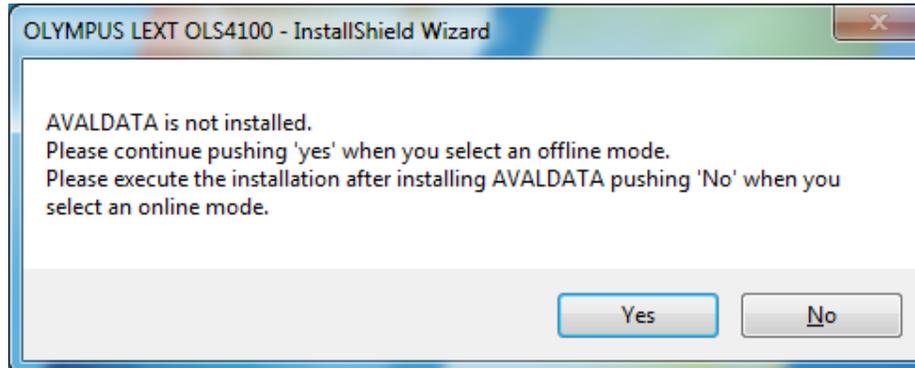


Fig.4 AVALDATA message

Select [Yes].

(This screen appears when the exclusive driver for I/F Board is not installed.)

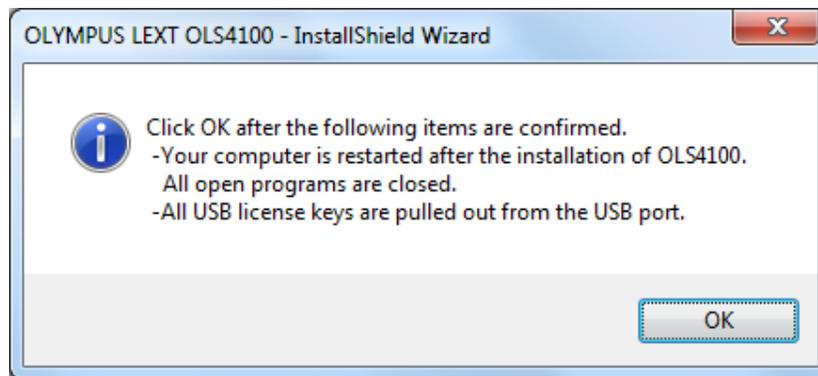


Fig.5 Confirmation message before the installation start.

Click [OK] after the following items are confirmed.

- As your computer is restarted after the installation of OLYMPUS LEXT OLS4100, all open programs are closed.
- All USB license keys are pulled out from the USB port.



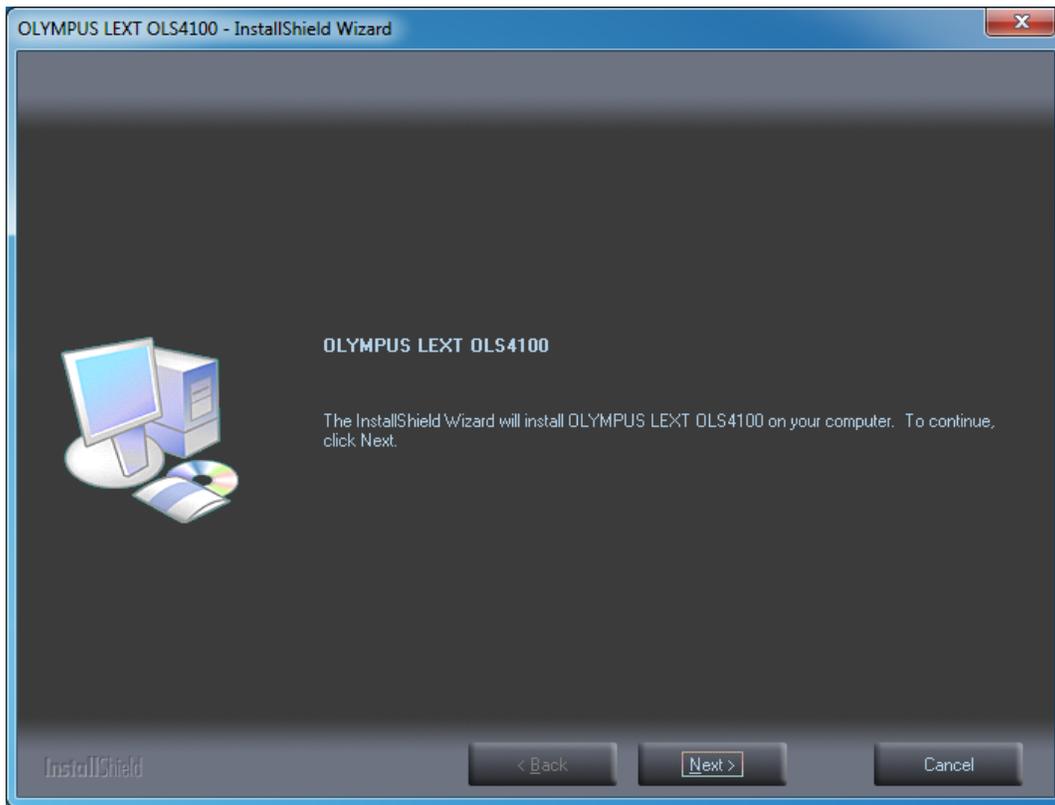


Fig.6 Welcome screen of OLYMPUS LEXT OLS4100

Select [Next].



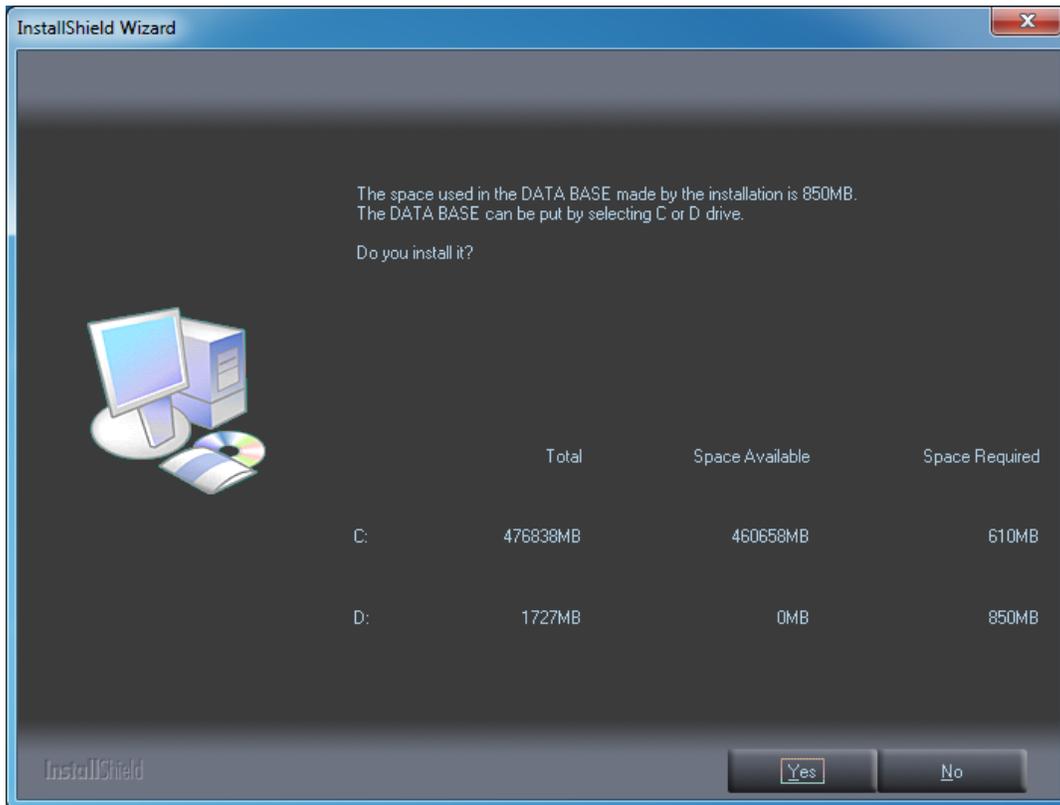


Fig.7 Confirmation of Hard Disk Capacity

Select [Yes]. Selecting [Yes] will start installing OLYMPUS LEXT OLS4100 Application (post-processing software).

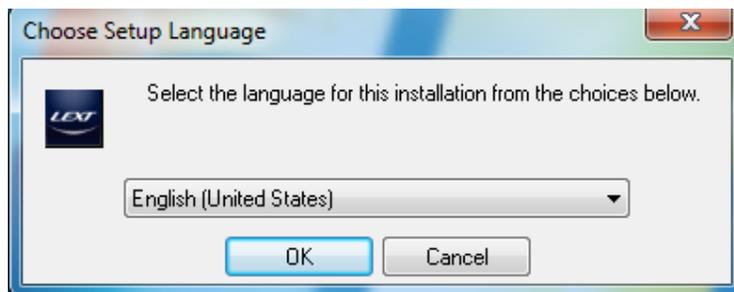


Fig.8 Selection of setup language in OLS4100 Application

Select English (United States), and select [OK].



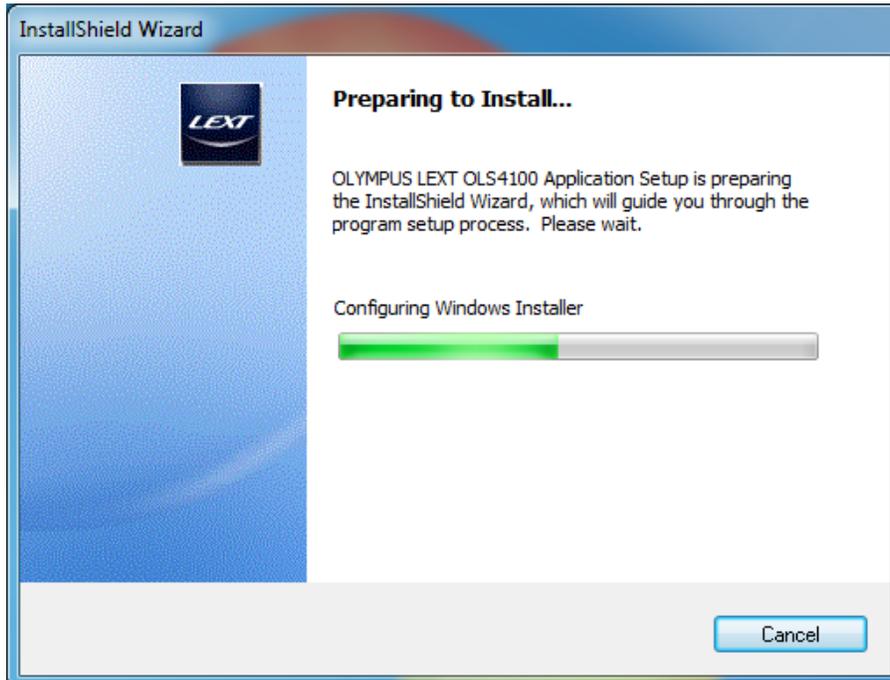


Fig.9 Preparation 1 to install OLS4100 Application

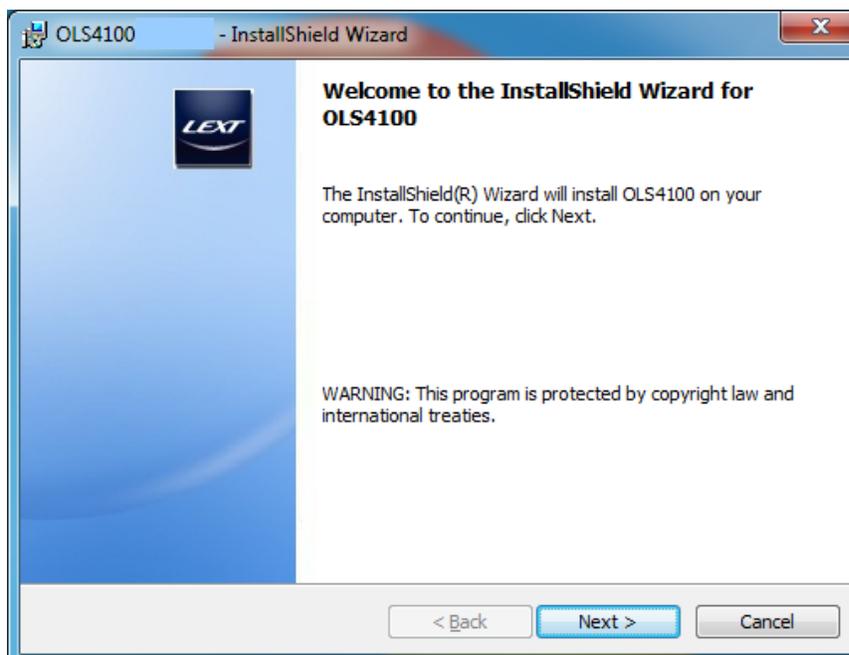


Fig.10 Welcome screen of OLYMPUS LEXT OLS4100
Select [Next].



Fig.11 License Agreement

Select **[I accept the terms in the license agreement. (A)]**, and select **[Next]**.

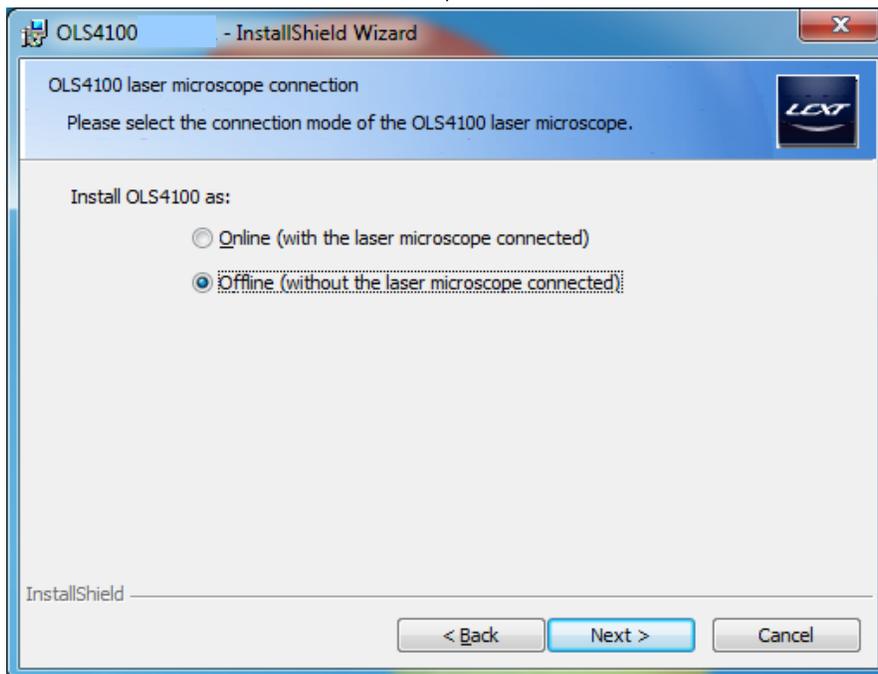


Fig.12 Connection Mode

Select [**Offline (without the laser microscope connected.)**], and select [**Next**].



OLS4100 administrator setup
Please enter the information about the administrator of OLS4100.

User ID
ADMIN

User name
Administrator

Password
olympus

Confirm password
olympus

InstallShield

< Back Next > Cancel

Fig.13 Creation of OLS4100 Administrator

Enter **olympus** in [Password] and [Confirm Password].

Select [**Next**].



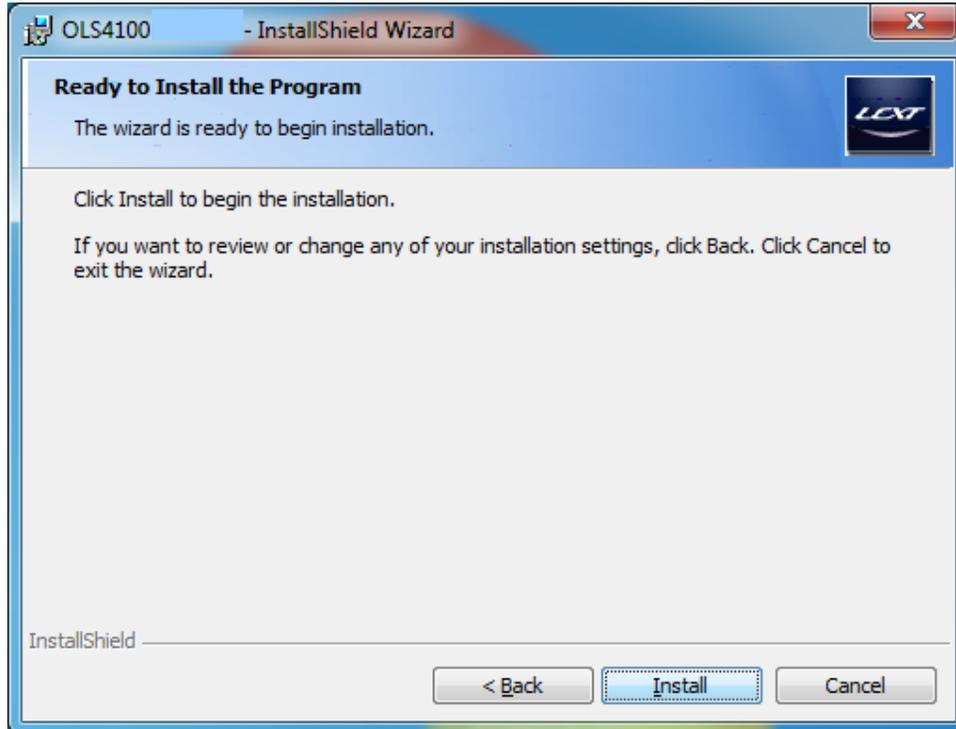


Fig.14 Preparation to install OLS4100 Application

Select [**I**nstall].



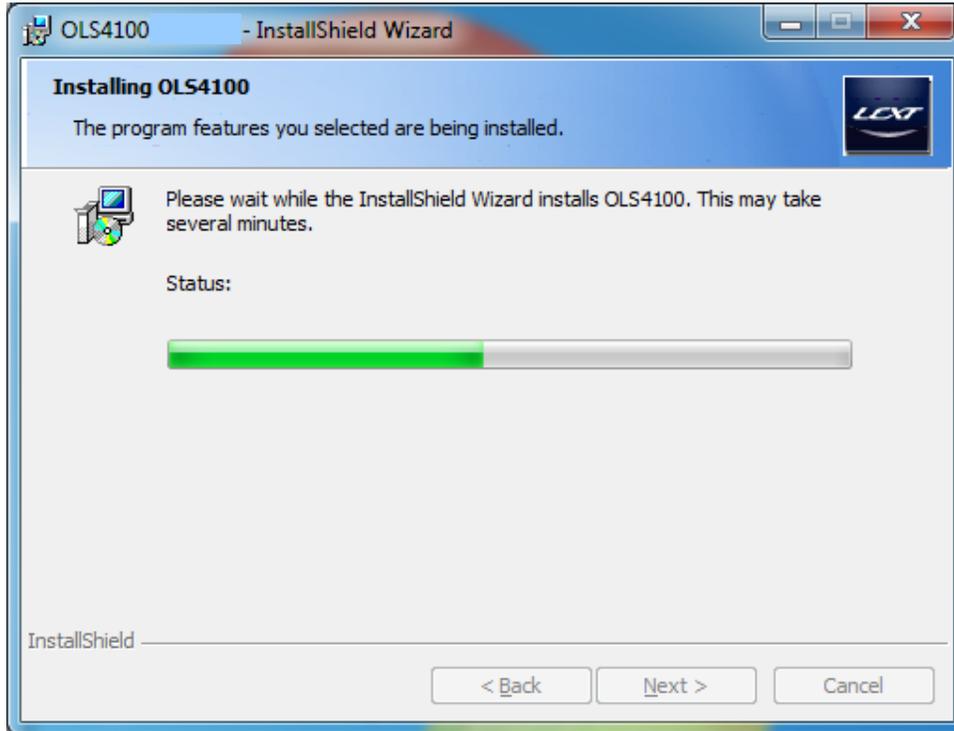


Fig.15 Installation of OLS4100 Application starts.

*** Please do not click "Cancel" button or "X" button of above installation window during installation. The controller environment state may not return to before installation start state.**



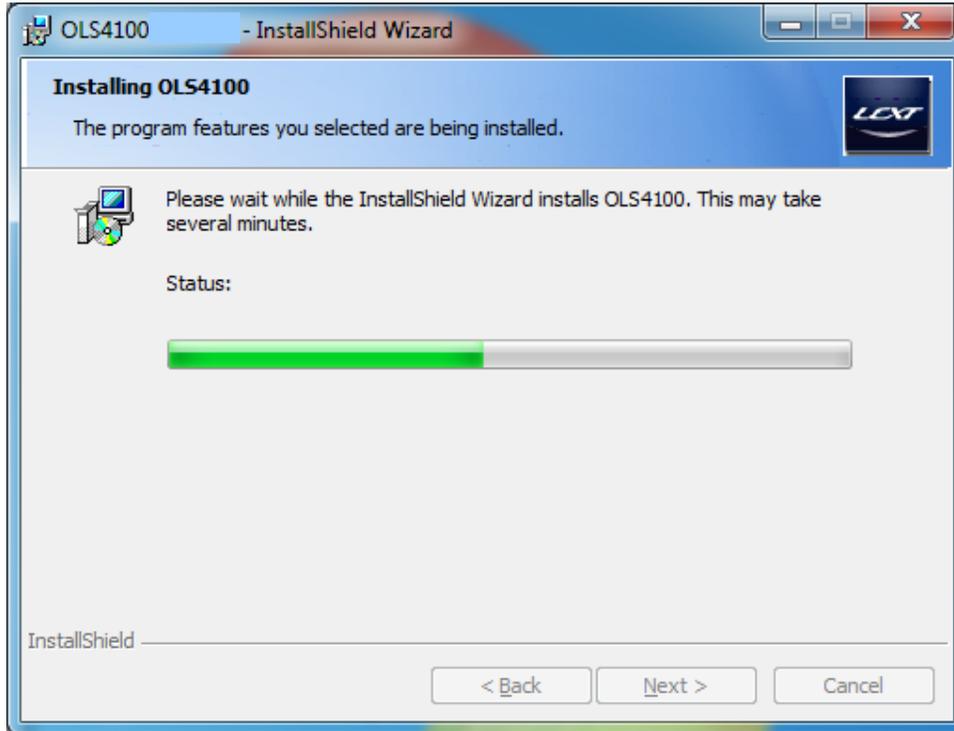


Fig.16 Installation of OLS4100 Application



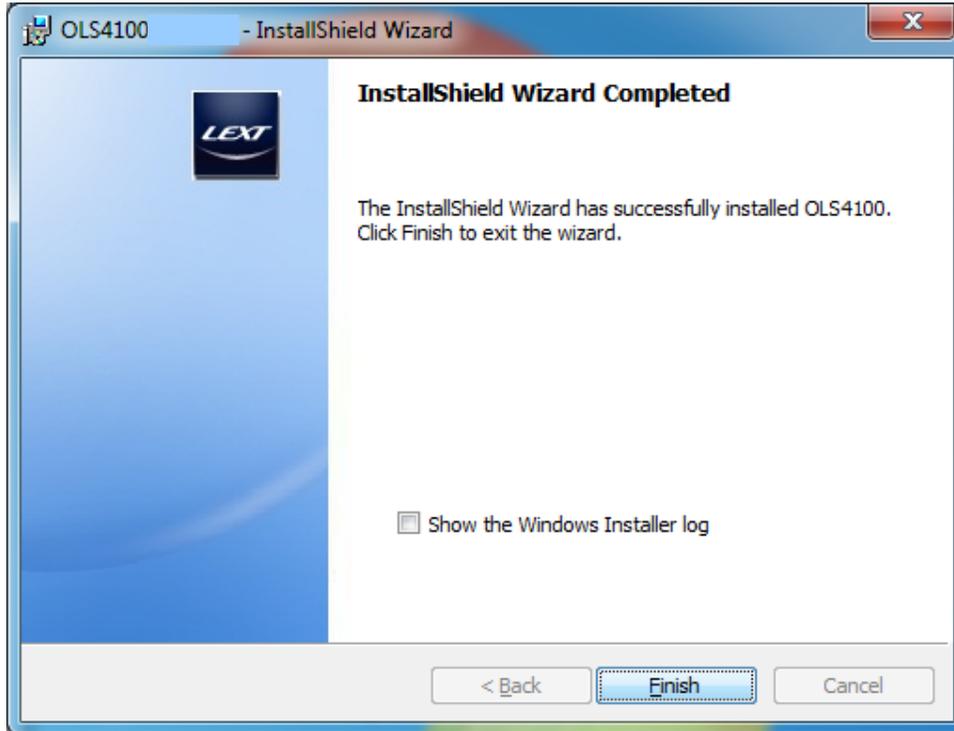


Fig.17 Installation of OLS4100 Application is successfully completed.

Select [**F**inish].

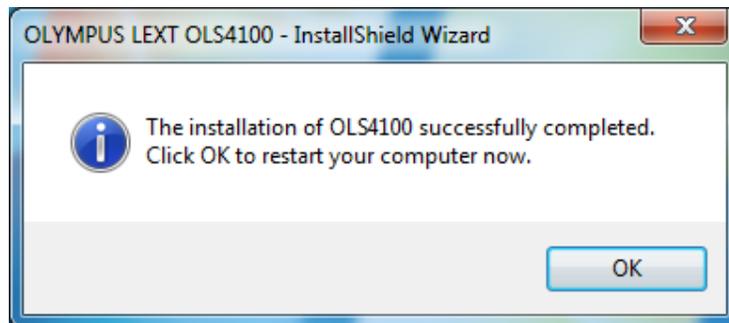


Fig.18 Completion of installation is displayed.

Select [**O**K] to restart the computer.

Warning Messages

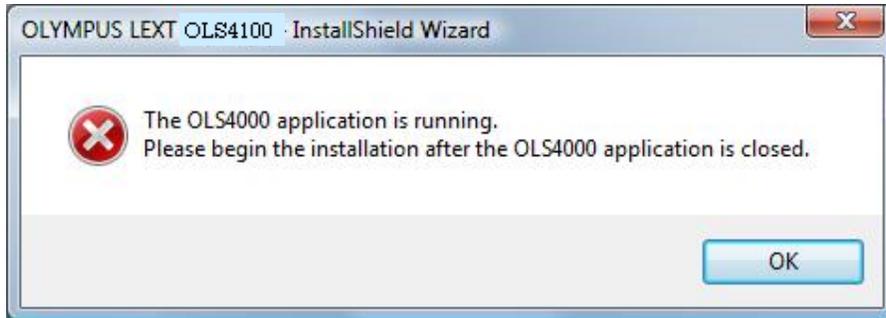


Fig.19 Warning for Application startup

This message appears if installation starts while OLS4100 Application is running. Close OLS4100 Application, and start installation again.

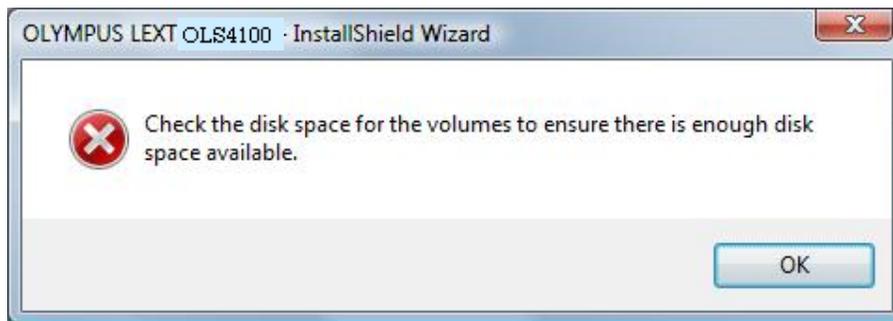


Fig.20 Insufficient Capacity

This message appears if the capacity of Hard Disk which is the destination of installation is insufficient. Secure the capacity, and start installation again.

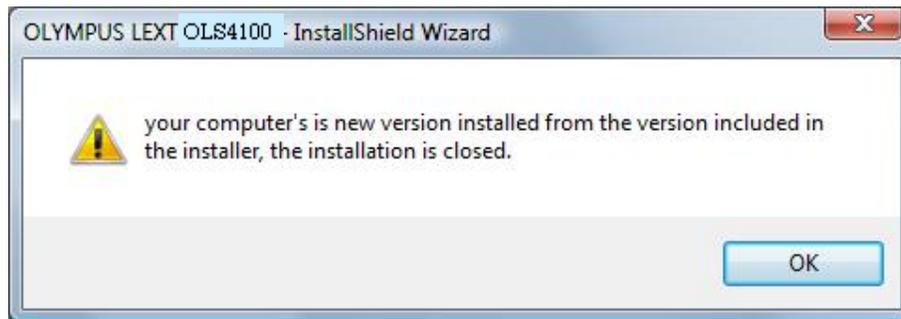


Fig.21 Display of Version check

This message appears if the version installed is newer than the version which is to be installed from now.

If you still want to install, uninstall the application installed now, and start installation.

End