
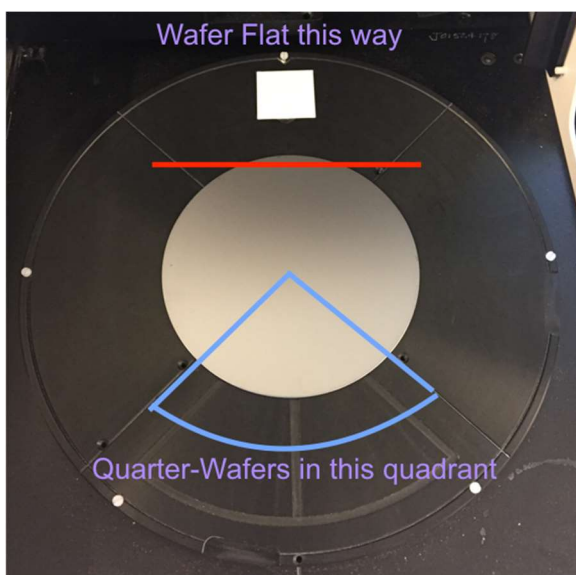


CDE ResMap Operating Instructions

Do not attempt to move your sample while the tool is running.

Probe conditioning may need to be performed periodically and is available under the “Operator” menu.

1. The tool is often left with the previous run's data open. If this is the case perform the following steps, if not proceed to step 5.
2. Select “Continue” on the upper left portion of the tool bar to exit the previous user's data screen.
3. Select “Continue” again.
4. Select “Cancel” on the lower middle portion of the software.
5. Load your sample onto the chuck in the manner illustrated below.

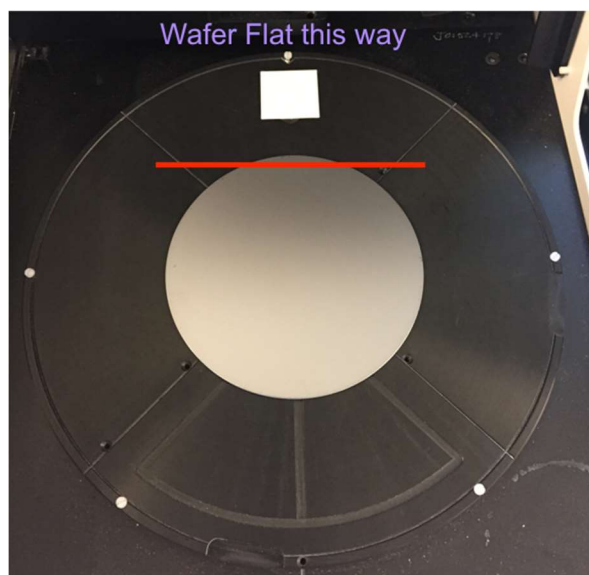

6. Select “Operator” from the tool bar on the upper left portion of the software.
7. Select “Run Recipe”
8. Select “UCSB_USER”
9. Select the appropriate recipe for your sample (in example, “4in_round 19pt”).
10. Select “Ok”.
11. Select “Run”.
12. After your run is complete, confirm your data.
13. Select “Continue” on the upper left tool bar to advance to the next data screen.
14. Select “Continue” again to leave the data display screen.
15. Select “Cancel” on the lower portion of the software to exit your recipe and return to the main menu.
16. Remove your sample.