

Substrate Template Name	Sub Size	Manually Find Center with Overview Camera
_Automatic small_OptAF*	< 12 mm	Yes
_Automatic rectangular_OptAF	>12 mm	No
_Automatic round_OptAF	>25 mm	No
<b>Separator</b>		
_Automatic small_PneuAF**	12 mm - 25 mm	Yes
_Automatic rectangular_PneuAF	> 25 mm	No
_Automatic round_PneuAF	> 25 mm	No

**\*RUNNING PIECES SMALLER THAN 5 MM REQUIRES ADDITIONAL TRAINING!!  
CONTACT STAFF**

**\*\*PIECES SMALLER THAN 12 MM CANNOT BE RAN WITH PNEUMATIC FOCUS!!  
SYSTEM DAMAGE WILL OCCUR!!**

**This is 12 x 12 mm →**

