






Blue M Oven Instructions





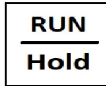


- Turn on: Middle switch below the door.

Other buttons:  The Back Arrow will cycle through the menu options: off, ctrl, PEnt, and P1-8.

- Cycle to desired recipe to use or edit (Example: P1), press the down arrow . This will load and start the P1 recipe.

- To stop the run for editing, immediately press the  Run/Hold button, then press the back arrow  button to cycle to “Off”, then press the down arrow .

Editing instructions:

1. Cycle in the menu to PEnt, then press the down arrow .
2. Then using the down arrow , cycle through the program’s options. To edit an option, press the back arrow  button.
3. Use the up and down arrows   to change an options value, and use the back arrow  to save the change made.
4. When done editing, use the up arrow  to get back to the main menu.

Program Options with typical values:

PN: 5
n5: 4
PLet: 1
dhru: 0
dhrd: 0
PEnd: 0

Specific program steps:

rt: Ramp Time
sp: Soak Point (temperature)
E1: on
E2: off
st: Soak Time
E1: on
E2: off